

Phage Lambda DNA EcoRI digest



0.7 % Agarose

bp	ng/1 µg
21226*	438
7421	153
5804	120
5643	116
4878	100
3530*	73

Features: Supplied in 6X Loading buffer

Description/Preparation:

The EcoRI digest of phage-DNA yields the 6 discrete fragments. Lambda DNA was completely digested by EcoRI, phenol extracted, ethanol precipitated, dissolved in 10 mM Tris-HCl (pH 8.0) and 1 mM EDTA. **The marker is ready-to-use.**

Usage: 0,5 µg/lane

Concentration: 0,4 µg/µl

Number of bands: 6 21226*, 7421, 5804, 5643, 4878, 3530*

Loading:

Agarose Gel / Polyacrylamide Gel

- Vortex gently before using
- apply 0,5 µg (agarose) or 0.5-0.9 µg (polyacrylamide gel) per 1 mm lane

Quantification:

See the graph for the percentage of the bands and the amount of DNA per band in ng, relating to 0.5 µg loaded marker. Use the same volume of DNA and marker. Additionally the concentration of loading buffer in samples and marker should be equal.

Note:

Dilute in TE or other buffer of minimal ionic strength. DNA may denature if diluted in dH₂O and subsequently heated.

* The cohesive ends (12b cos site of bacteriophage) of fragments 21226 bp and 3530 bp may anneal to and form the additional band. These fragments may be separated by heating to 65°C for 5 min and then cooling on ice for 3 min.

Transportation: Shipped on blue ice or room temperature

Storage: at -20°C for 24 months

Ordering information:

Cat.-no	Description	Amount
300011	Phage Lambda DNA EcoRI	2 x 100 µg
300012	Phage Lambda DNA EcoRI	5 x 100 µg